

STK[®]

Sperm Tracker

SPRAY


ON THE RAPE SCENE
FOR HARD SURFACES

ISO 18385
Forensic Grade

STK SPRAY, DIRECTLY SCREEN THE RAPE SCENE

STK Spray is a must-have in CSIs' toolbox to screen directly the rape scene. If today clothes and bed sheets are often collected and sent to the lab, the rape scene is often not or poorly being screened due to poor results given by ALS for semen screening. Key traces are surely missed. How to be sure no evidences are missed? How to screen non-transportable evidences? How to save time and improve investigation success rate?


STK Spray is the first screening solution that enables a **sharp and specific localization of sperm traces directly on the rape scene** (floor, furniture, door handle, sink, tree leaves, grass, gravels...). STK Spray disrupts current procedures by **instantly** and **reliably** detecting semen traces directly on the scene. STK Spray is **ready-to-use**, just dissolve one pouch in water and you are ready!

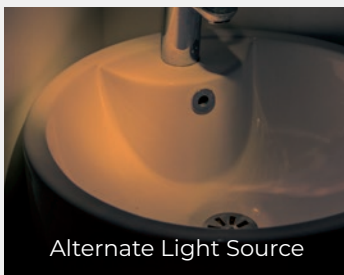

Alternate Light Source



STK Spray

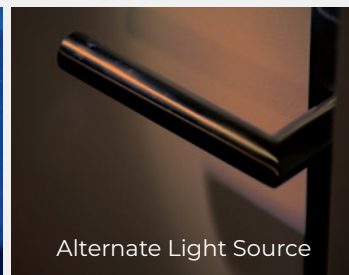

Alternate Light Source



STK Spray

STK Spray offers many advantages to CSI for semen stains detection


- **Sensitivity:** even mixed, wiped and rinsed traces will appear
- **Specificity:** only semen stains will be visible
- **Non Toxic:** human and environment-friendly
- **Do not alter DNA,** fully compatible with extraction techniques and DNA analysis
- **Convenient** and **portable** solution
- **Easy** and **ready-to-use:** dilute 1x STK pouch in demineralized water
- **ISO 18385** Forensic Grade & **ISO 13485** standards
- Box of **sealed individual pouches**
- 1 pouch ~10m²

EASY & RAPID PROTOCOL ON THE CRIME SCENE



1

Dilute 1 pouch of STK Spray
in 100 ml (~3 fl oz)
of demineralized water

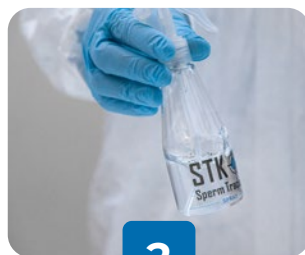


2

Shake lightly,
wait 30 seconds.
Solution is ready-to-use



3

Spray the scene
at a distance of ~20 cm
(~8 inches)

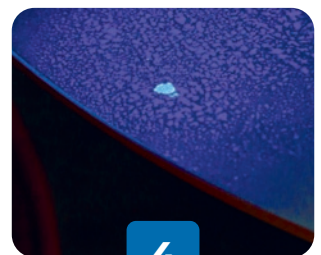

4

Read results with STK UV
Light: only semen stains
will glow in a blue color



STK SPRAY

- Can be used indoor and outdoor
- **100% of semen stains are detected** on the field
- **No false positive** and above all **no false negative**

AXO-STK-SP-10: STK Spray - Box of 10 pouches

SPRAY BOTTLE

- Fine mist
- Perfect to spray STK Spray

AXO-STK-PV1

POSITIVE CONTROL

- Box of 10 positive controls
- Perfect to validate your STK Spray solution

AXO-STK-PC-10



STARTER KIT STK SPRAY

- STK Spray: box of 10 pouches
- PC10: box of 10 positive controls
- Rechargeable STK UV Light
- Spray bottle
- UV glasses (clear)
- Case

AXO-STK-SPRAY-STARTER

STK UV LIGHT



- Portable light with battery.
Rechargeable, perfect to bring on the field

AXO-STK-L-UV

AXO Science

AXO Science develops innovative solutions in the field of molecular biology and forensic. We are now bringing to forensic scientists the STK Sperm Tracker technology, which is revolutionizing the detection of sperm traces on the rape scene and in the laboratory. AXO Science is ISO 13485 and ISO 18385 certified.

Contact us

sales@axoscience.com
www.sperm-tracker.com
Follow us on  

For forensic use only

Copyright © 05/2025 AXO Science SAS. Non contractual document. All rights reserved.

AXO Science, STK Sperm Tracker and related logos are trademarks or registered trademarks of AXO Science SAS. Manufactured by AXO Science SAS. (Ref) 25/05/AXOSTK/SPRAY/PM/0001-EN